




These installation instructions describe how to upgrade the RoboSep™-S Software Version 1.0.4.0 to RoboSep™-S Software Version 1.1.0.4. A keyboard, mouse, and USB key are required for this task. Please note that software version 1.1.0.4 is only compatible with the RoboSep™-S instrument. Do not attempt to install this update onto a RoboSep™ (Classic) instrument.


- 1 Visit www.stemcell.com/instruments to download the RoboSep™-S software. Save the file onto a USB key.
- 2 Plug in a keyboard and mouse into RoboSep™-S. Insert the USB key with the saved file into RoboSep™-S.
- 3 On the RoboSep™-S, go to the 'Maintenance' screen and press the 'Exit to Desktop' button. Press the 'Window' key on the keyboard, access 'My Computer' folder and select the USB key. Locate the saved file on the USB key and copy the file onto RoboSep™-S.
- 4 Run 'Update_RoboSep-S.exe'.
- 5 Select 'Yes'.

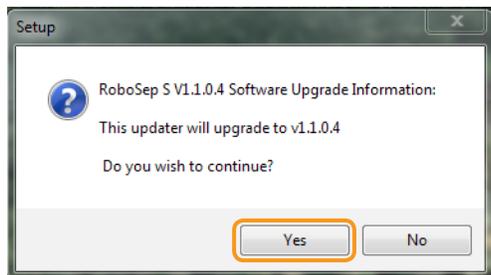


- 6 When the following dialog box is shown, click 'Next'.




- 7 Then click 'Install'.

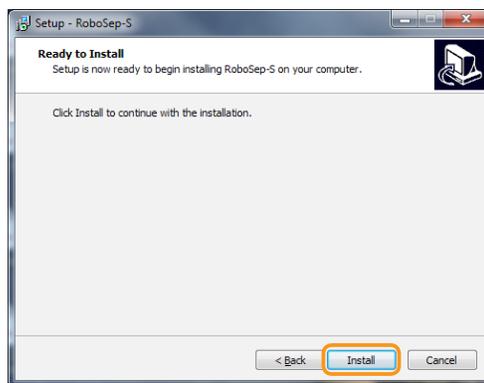

- 8 To upgrade the RoboSep™-S Server, select 'Next'.

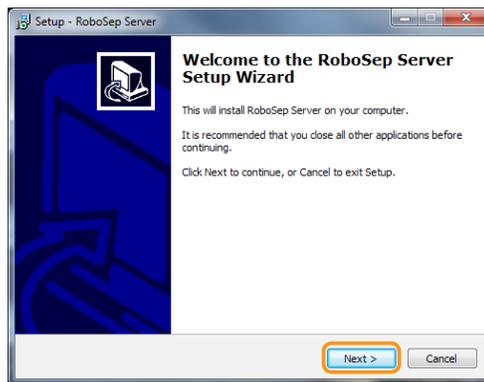


- 9 Continue by selecting 'Next'.

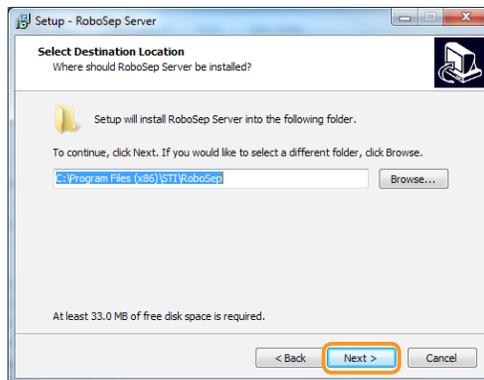




- 10 At this screen, select 'Manual Configuration' and then click 'Next'.

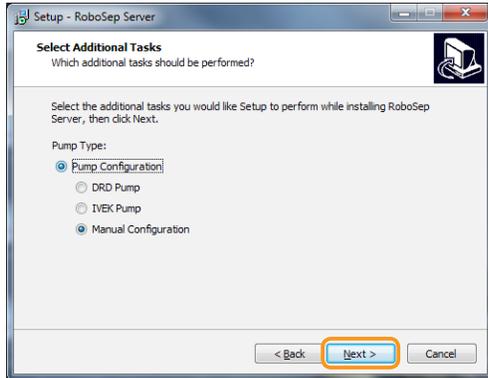


- 13 RoboSep™-S Server upgrade is now complete. Select 'Finish'.




- 11 Click 'Install' to start the installation.

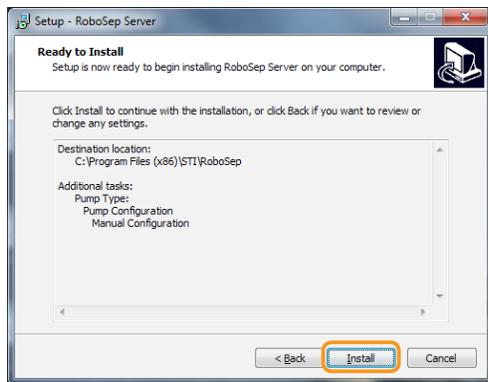


- 14 Continue by selecting 'Next'.

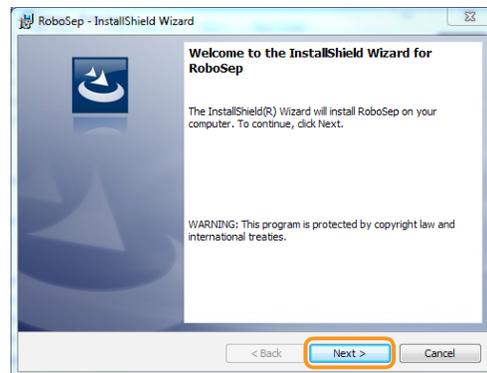

- 12 When this dialog box pops up, select 'OK'.

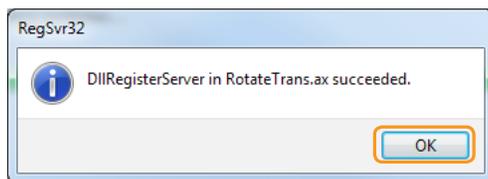


- 15 Continue by selecting 'Install'.

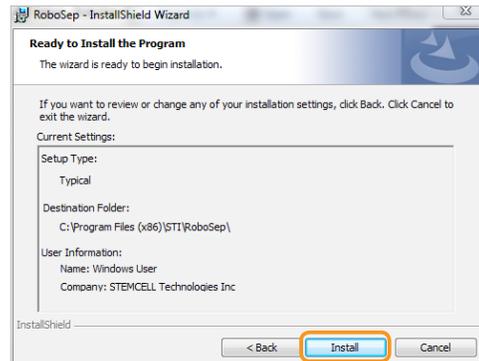





16 Click 'Finish'.

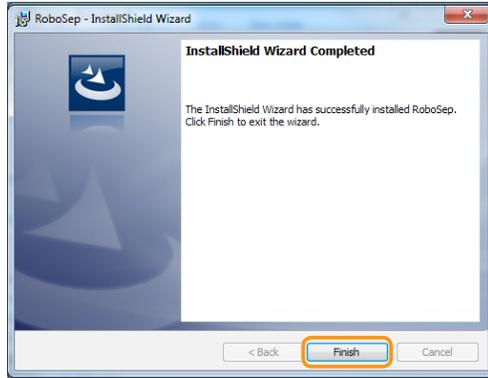


17 To upgrade the RoboSep™-S Service program, select 'Next'.

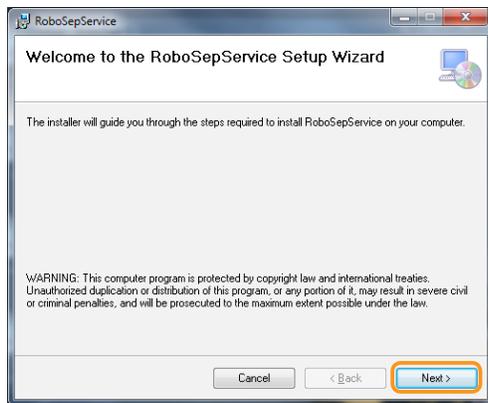


18 Upgrading RoboSep™-S Editor is started. Click 'Next'.

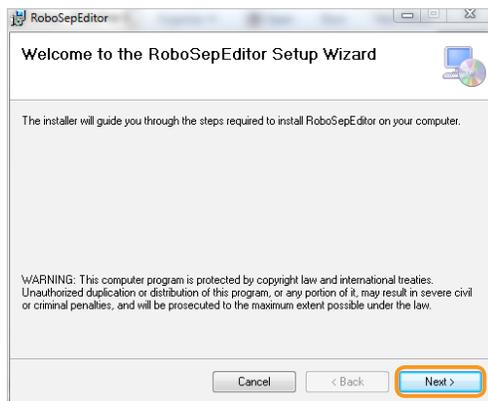


19 Click 'Next' when you see the following dialog box.

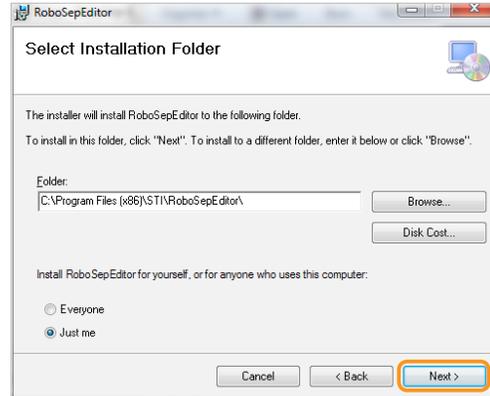

20 Click 'Next'. Copying files will start.

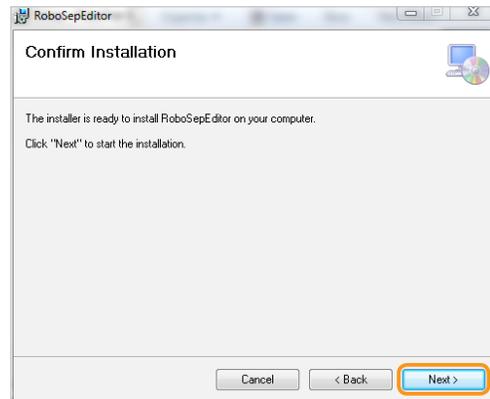

21 Upgrading RoboSep™-S Client is completed. Click 'Close'.

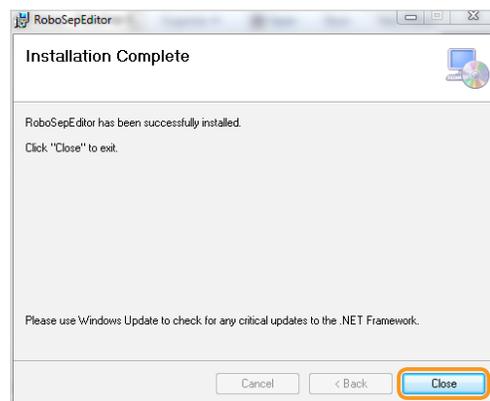





- 22 Select 'Next' at the following screen.




- 23 Select 'Next' at the following screen.

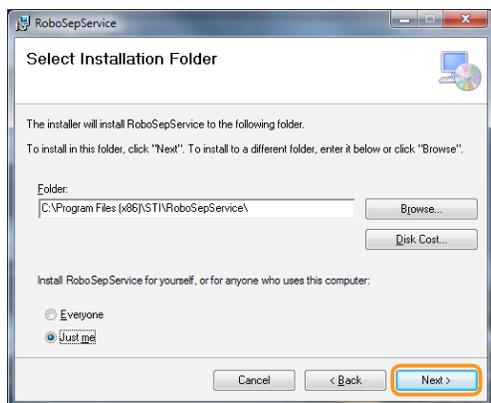


- 24 Click 'Next'. Copying files will start.

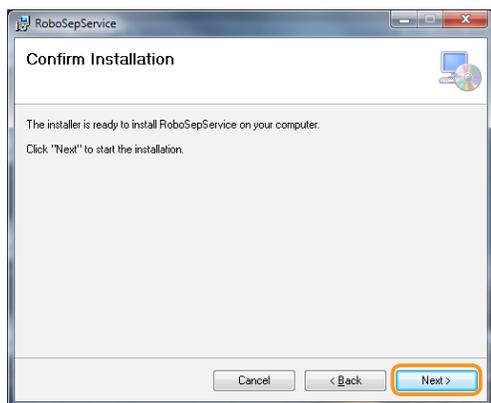

- 25 Click 'Close' at the following screen. Upgrading RoboSep™-S Service is completed.

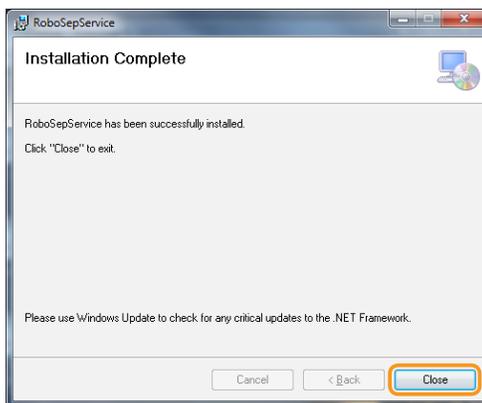

- 26 Click 'Finish'.

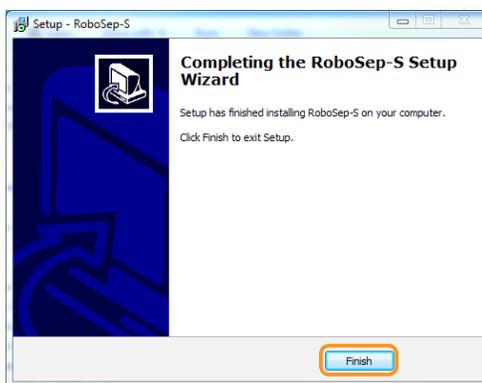

- 27 Press the Window key on the keyboard to restart your RoboSep™-S and return back to the RoboSep™-S Run Samples screen.