


Install Instructions for STEMvision™ Software Version 2016.04.21.00

These install instructions describe how to upgrade STEMvision™ Software Version 2015.12.18.01 to STEMvision™ Software Version 2016.04.21.00. An Internet connection, USB key, keyboard and mouse are required for this task.


- 1 Visit www.stemcell.com/instruments to download the newest STEMvision™ software. Save the file onto a USB key.
- 2 Insert the USB key with the saved file into the computer connected to the STEMvision™ instrument.
- 3 Run 'STEMvisionUpdate.exe'. Ensure dongle is inserted.
- 4 Select Next.




Note: This update is for the STEMvision™ Color Unit (Catalog #22000C). If you attempt to perform this install on a computer that is not connected to STEMvision or is connected to another unit, the following error is shown. Please contact your STEMCELL representative if you receive this error message.

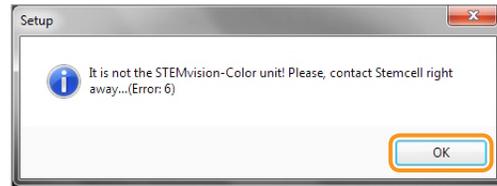


- 6 Ensure C:\CFC is entered as the installation path, and select Next.

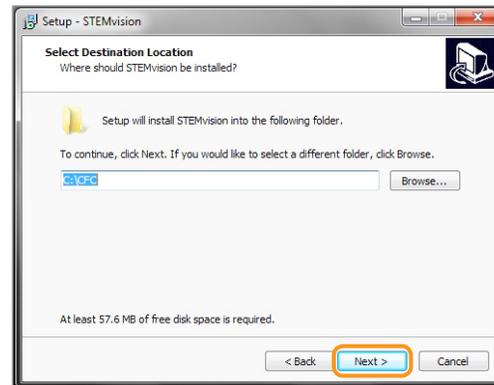


- 5 Enter the serial number of your instrument and select Next.

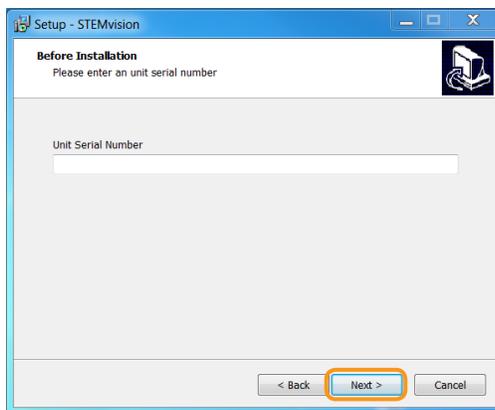

- 7 Continue by selecting Next.

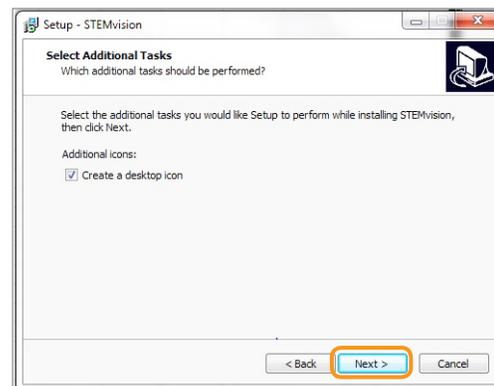

8 Select Install.

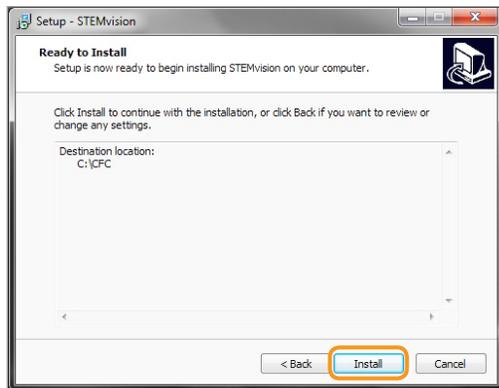

10 The STEMvision™ update is now complete. Select Finish.



9 Wait while the installation occurs.

