



Release Notes

STEMvision™ Software Version 2021.05.07.00

These release notes describe the new functionality and/or fixes to the previous STEMvision™ Software Versions 2018.04.16.00

Software	New Functionality/Fix
STEMvision™ Acquisition Version 2.5.0.1	<p>Data security enhancements:</p> <ul style="list-style-type: none"> Improved event logging of instrument application usage Added password protection to camera manual settings Option for data authenticity check Option for encryption of application log files <p>Bug fixes for improved stability.</p>
STEMvision™ Analyser (Human) Version 2.5.0.1	<p>Data security enhancements:</p> <ul style="list-style-type: none"> Improved event logging of instrument application usage Database password protection Option to add user level restrictions for analysis and report generation Option for data authenticity check Option to encrypt application log files Option for e-signature on PDF reports <p>Bug fixes for improved stability.</p>
STEMvision™ Analyser (Mouse) Version 2.5.0.0	No updates
STEMvision™ Colony Marker Version 2.5.0.1	<p>Data security enhancements:</p> <ul style="list-style-type: none"> Improved event logging of instrument application usage Option for data authenticity check <p>Added capability to comment on individual colonies</p> <p>Scale bar on image preview</p>

Copyright © 2021 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and STEMvision are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO WWW.STEMCELL.COM/COMPLIANCE.

Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713 • INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM
 FOR GLOBAL CONTACT DETAILS VISIT WWW.STEMCELL.COM DOCUMENT #27197 VERSION 1.0.0 JUN 2021